

Item Name : Cartesian Diver

Item Code : 150.2008.01

Item Description :

Cartesian Diver is supplied with hollow glass diver, with a 10" x 1.5" glass cylinder, a rubber sheet and rubber band. Fill the cylinder with water, drop in the diver so he floats, fasten the sheet totally across the top and then gently depress the sheet. The diver will dive, his speed depending upon the pressure you place on the sheet, release the sheet and the diver returns to the surface. The diver's behaviour illustrates the uniform transmission of pressure in all directions, the compressibility of gases and Archimedes Principle.

Item Image :



Cat. No.

150.2008.01